Inversina Taumelmischer 50 L/100 L/350 L

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Inversina tumbler mixer 50 L/100 L/350 L

The Biocomponents inversina tumbler mixer uses inversion kinematics for mixing of solids and liquids. This astonishing three-dimensional motion (Paul Schatz principle) combines a figure-of-eight movement with rotating and sets the substances in a rhythmic, pulsating motion.

Characteristics

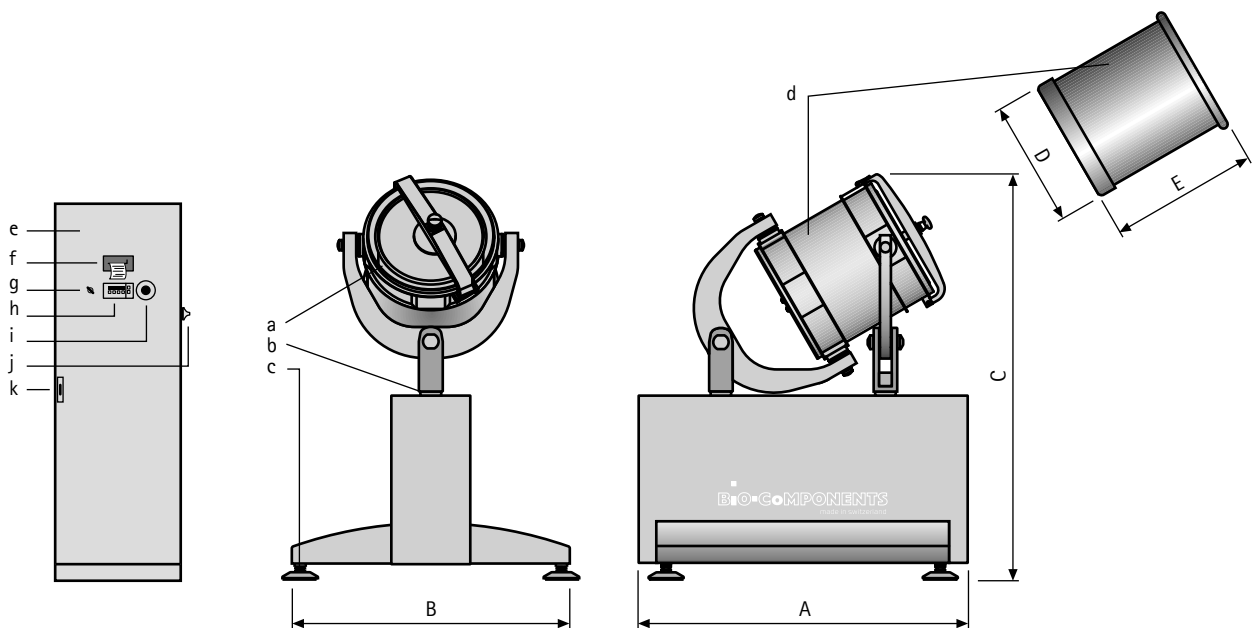
- Intensive, fast and very gentle mixing of components of different density, size, shape and concentration.
- Best mixing quality and no segregation due to the inversion kinematics.
- High quality mixing results independent of vessel shape and without internal agitators.
- Dust-free and hygienic mixing in closed containers.
- Two driven axles.
- Easy cleaning and maintenance.
- On request adapted to customer specific containers.
- Safety concept and mixing equipment as independent units.
- Control unit located outside of safety zone.
- Available also in GMP design.

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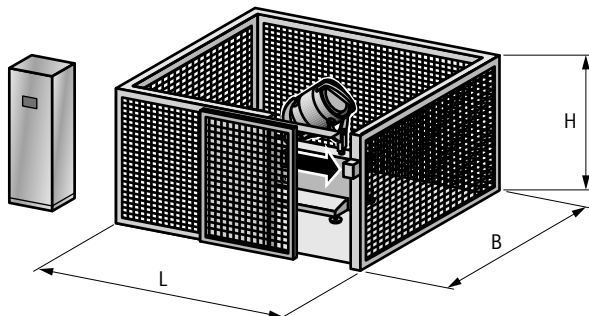
Safety concept

To assure the protection of personnel, the Inversina must be operated within a safety zone. Bioengineering offers an optional device with a safety door designed for permanent installation.

The Inversina can only be started if the safety door is closed. During operation, as long as the Inversina is in motion, the safety door cannot be opened. The latch is not released until the Inversina has stopped; then, the safety door can be opened.

For unhindered operation, the control unit of the Inversina is set up outside the safety curtain.

If the structural situation or other circumstances prevent the installation of such a safety curtain or if the Inversina is to be operated in a separate room, the buyer shall be responsible for the safety concept and wiring. Bioengineering cannot accept any liability for solutions that deviate from the standard safety configuration.

L x B x H [mm]

50L: 1345 x 1345 x 2116

100L: 1680 x 1680 x 2030

350L: 2470 x 2070 x 2516

L x B x H [inch]

50L: 52.95 x 52.95 x 83.31

100L: 66.14 x 66.14 x 79.92

350L: 97.24 x 81.50 x 99.05


Safety curtain

Order no.

Safety curtain, 50 L, Plexiglass	*
Safety curtain, 50 L, Steel grid	*
Safety curtain, 50 L, Steel sheet	111221
Safety curtain, 100 L, Plexiglass	13159
Safety curtain, 100 L, Steel grid	*
Safety curtain, 100 L, Steel sheet	*
Safety curtain, 350 L, Plexiglass	13271
Safety curtain, 350 L, Steel grid	13602
Safety curtain, 350 L, Steel sheet	13568

* on request

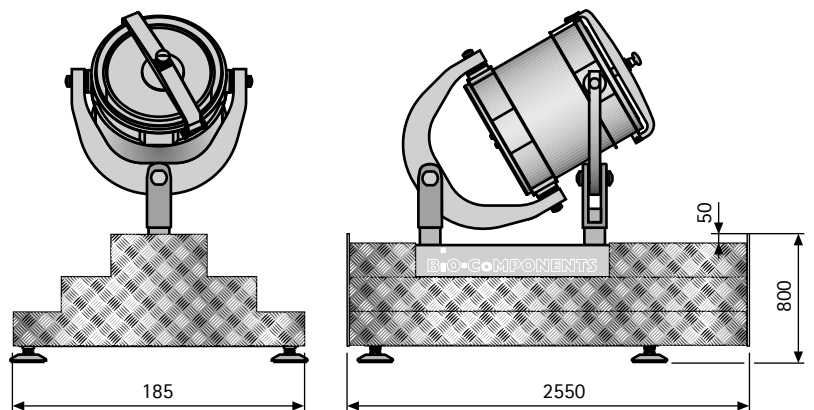Operator stair

Order no.

Operator stair	51416
----------------	-------

Steel container (Material 316 L)

Order no.

50 L	52040
100 L	52041
350 L	52044

Spare parts

on request